

Install for windows7 32bit&64bit

Note that 32bit and 64bit are different driver software

You should confirm your PC as which bit type.

Insert the “Mpixel Driver for windows 7 “ CD to your PC.

Caution : Do not connect USB cable to your PC

Open the “IUC Driver” in the CD

Select 32bit or 64bit

Appear as below




Windows 7 32bit




Windows 7 64bit

Double click the “dpinst “, and automatically

32ビット

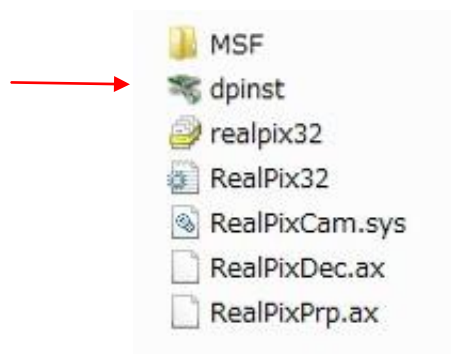

64ビット

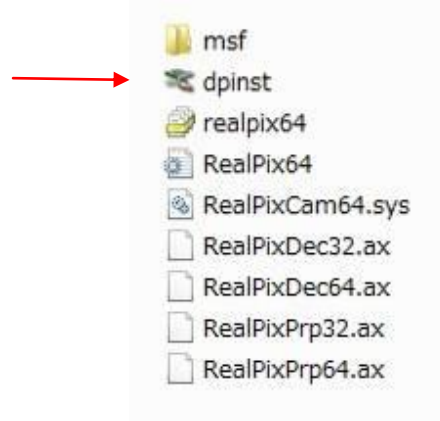


Start as below image


Click the “NEXT”



Starting the install

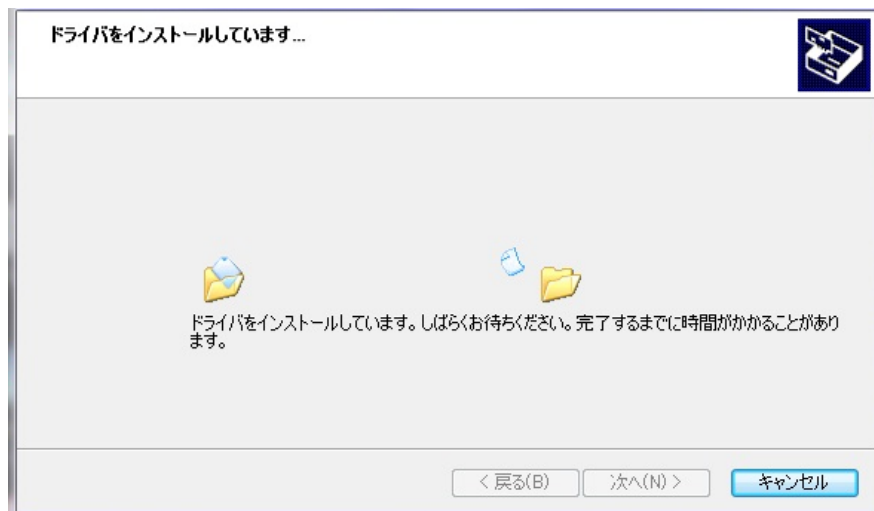

Click "Install"

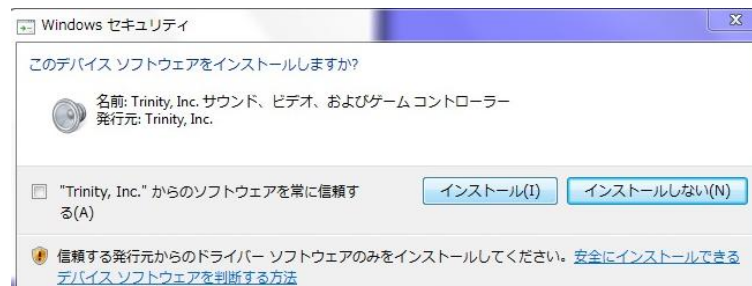

Click "Finish"

。



After finished the install, you can connect the USB cable to your PC.